



- Remove the rivet on the door hinge side in the pull-out shelf by drilling. (Fig. 3)
- Position the locker.
- Align the holes on the door hinge side with each other.
- Connect locker and pull-out shelf with rivet supplied. (Fig. 4)

- Warning!**
- The self-closing of the safety storage cabinet must not be impaired by the installation of a locker:
- Install the locker so that the door hinges of the locker door and the cabinet door are on the same side. (Fig. 1)
 - For double-leaf doors, install the locker on the left or right side. (Fig. 2)

Parts list		
Pos.	Title	Quantity
010	Locker	1
020	Rivet	1

Article number Locker	29-50000-037(L)
-----------------------	-----------------

This drawing is our property! Furnish this information to other persons and reproduce or copy is not permitted without our written permission.	title: Installation instruction Locker CLASSIC pro / PREMIUM pro	
	drawing no.: 000123591	page: 1 of 1
	revision: 01	date: 2023-06-13
Düperthal Sicherheitstechnik Frankenstrasse 3 D-63791 Karlstein		