
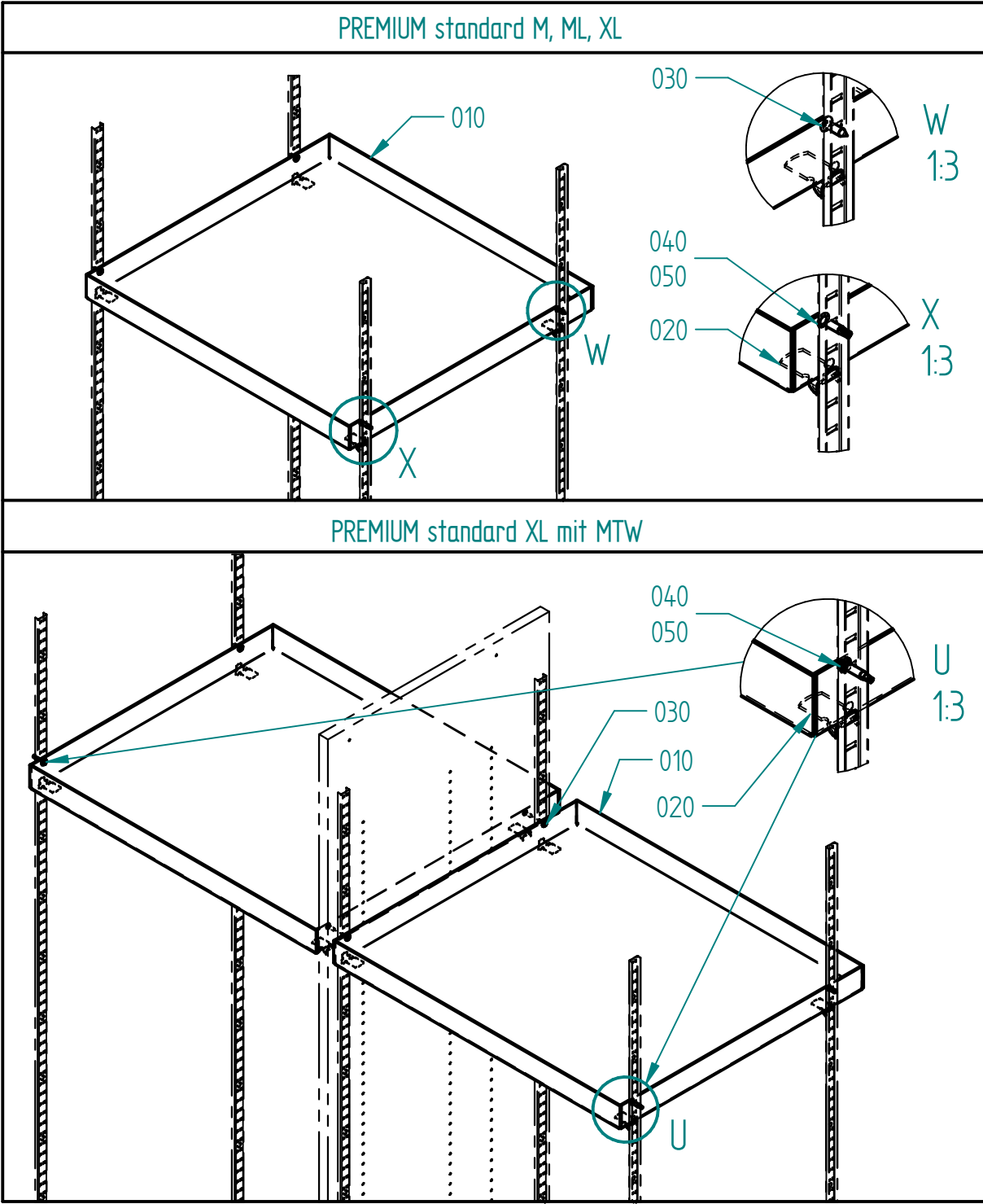


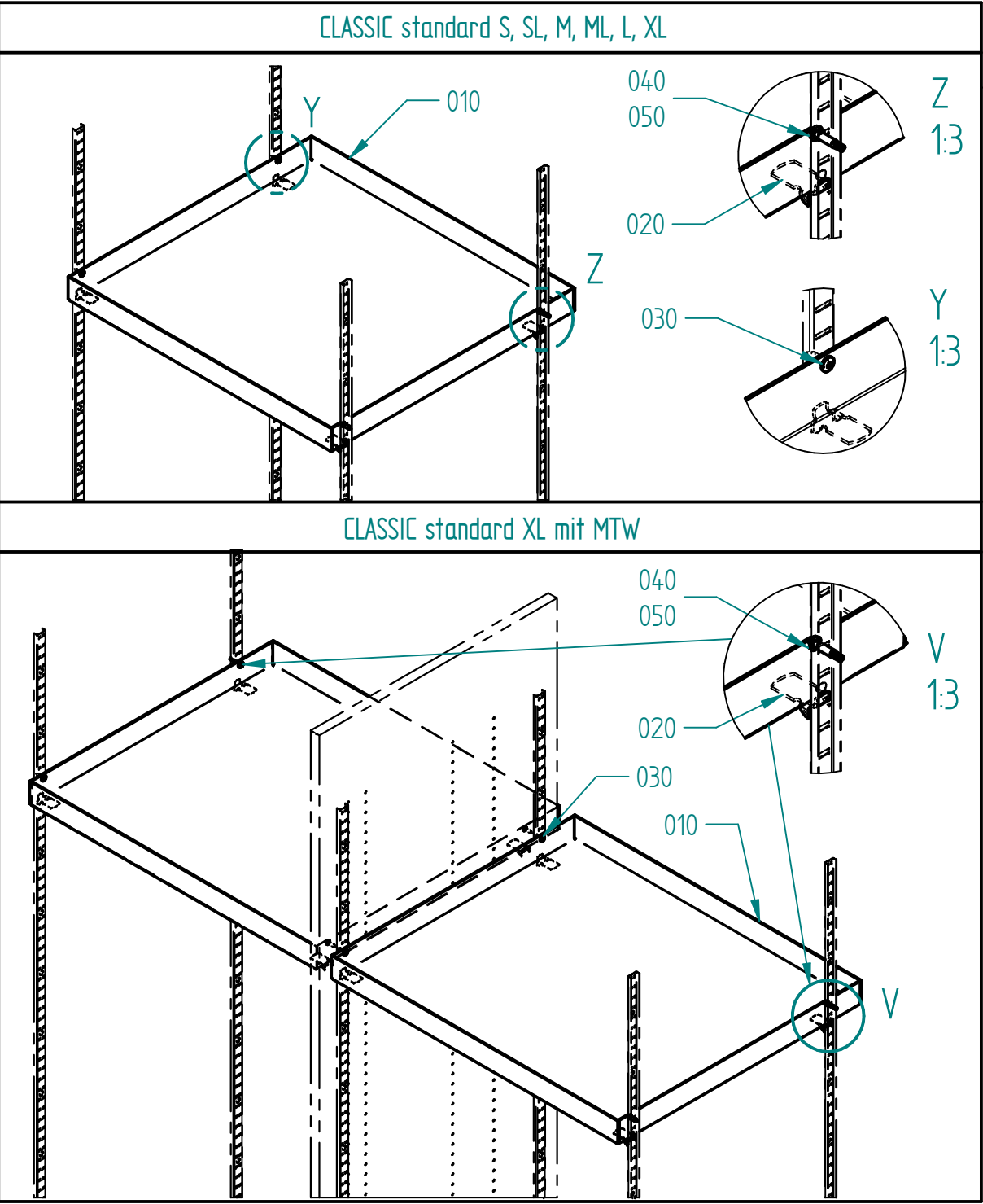

| Parts list |                           |       |
|------------|---------------------------|-------|
| Pos.       | Description               | Piece |
| 010        | storage shelf             | 1     |
| 020        | shelf support             | 4     |
| 030        | crosshead screw           | 3     |
| 040        | drill screw               | 1     |
| 050        | serrated lock washer      | 1     |
| 060        | installation instructions | 1     |

Installation of the additional storage shelf:

1. Insert shelf support pos. 020 into the shelf support rails at the desired height
2. Place storage shelf pos. 010 on the shelf supports
3. Screw the storage shelf to the shelf support rails using pos. 030, pos. 040 and pos. 050  
(Screw connection varies depending on the cabinet, see illustration)

This drawing is our property!  
Furnish this information to other  
persons and reproduce or copy is  
not permitted without our  
written permission.

Düperthal Sicherheitstechnik  
Frankenstrasse 3  
D-63791 Karlstein

title: Installation instructions  
for additional storage shelf

drawing no.: 000093859

page: 1 of 1

revision: 01

date: 2018-08-01

 **DÜPERTHAL**